

U-react Powder for Uranium Test

Each U-react vial contains enough U-react reagent to perform 150 tests. This Uranium test kit works for 5 mL samples.

Storage

Store between +2 °C and +25 °C, away from light. Expiration date on box label refers to unopened material. Once reagent vial is opened, use within six months.

Instructions

1. Refer to your Uranalyze® test kit instruction manual for sample preparation. At the point where the manual directs you to add reagent, proceed as outlined below.

- 2 Add a dose (0.1 g) of U-react powder into the sample (5 mL) with the measuring spoon.
- 3 Cover the sample container and shake to dissolve the U-react powder.
- 4 Proceed as described onto Uranalyze® instructions.
- 5 Dip the test strip for 60 seconds and take on immediate reading.
- 6 Compare the test field with the colorimetric scale indicated on the strip container.

